


User Manual
FlexiVap-24
Automatic Intelligent
Parallel Concentrator

More Intelligent, More Efficient

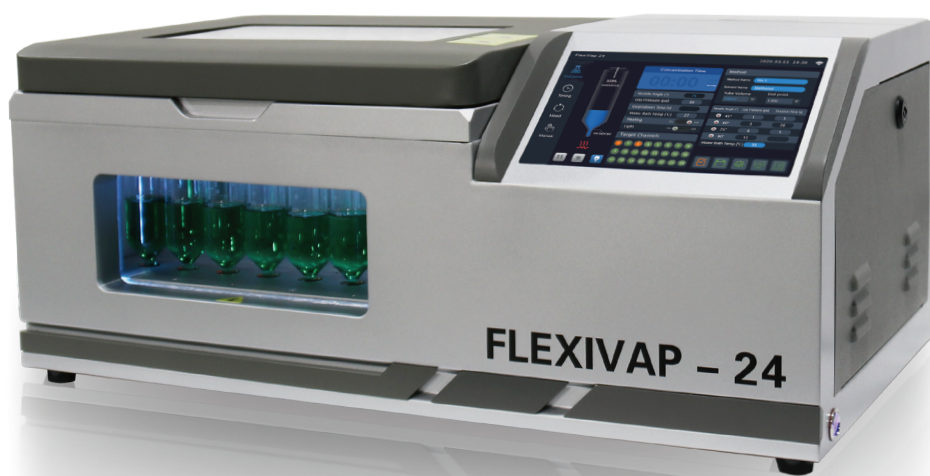

Cautions

FlexiVap-24 Automatic Intelligent Parallel Concentrator enables to adjust automatically the nozzle angle. Please do not pull the nozzles or the rotation axis manually, otherwise the equipment may be damaged.

Nozzles will automatically reset after powering on. During the concentration process, gas will stop blowing when the cover is opened, and start again with the cover closed.

Please be acquainted with the following icons before using this equipment.



Warning Danger (please be careful and note safety)



Key Information (please kindly notice)



Nozzle angle will be automatically adjusted, please do not operate manually.

CONTENTS ▶▶▶

Cautions	01
Contents	02
Announcements	03
I. Introduction	04
II. Unpacking Instructions	05
III. Installation	05
IV. Environment	05
V. Hardware	06
1. Control Terminal	06
2. Top Cover	06
3. Water Bath & Sample Rack	07
4. Exhaust Device	07
5. Power Socket & Button	07
6. Gas Inlet & Drainage Outlet	07
7. Lighting	07
VI. Software	08
1. Main Interface	08
1.1. End-point Mode	08
1.2. Timing Mode	08
1.3. Mixed Mode	09
1.4. Manual Mode	09
2. Method Settings	12
VII. Operation	13
1. Method Settings	13
2. Loading & Application	14
3. Process Pausing	14
VIII. FAQ & Solutions	14
IX. Service	15
Statement	16